Manufactured By: HYPHEN BioMed

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CONSERVER A  
2-8 °C

**Origin:**

Highly purified factor X from citrated bovine plasma, activated with agarose bound RVV.

**Presentation:**

Vial containing approximately 15 µg of highly purified bovine factor Xa in presence of additives and preservatives.

(1 Unit FXa is the amount which can be generated from 1 ml citrated bovine plasma, ie about 10 µg).

**Reconstitution:**

Each vial must be restored with 6 ml distilled water, in order to get a buffered solution (pH7.90) at a FXa concentration of about 2.5µg/ml. It can be diluted to the desired concentration, when required, in the adequate buffer.

**Excipients:**

Content per vial: Tris , Sodium chlorure, PEG 6000, Bovine Serum Albumin, stabilizers.

**Purity:**

The factor X zymogen used for preparing factor Xa has one major band of about 55,000 daltons on SDS-PAGE.

**Activity:**

- a) Tested for its amidolytic activity on a FXa specific substrate, in a purified system without or with RVV. All FX is converted to FXa.
  - b) Tested in the anti-Xa heparin assay with substrate Xa.
  - c) Chromogenic activity on specific substrates (expressed in nkats):
- When tested in the optimized conditions (0.05M Tris buffer at pH8.40, containing 0.30M NaCl), the following specific activities are observed on the various FXa substrates (expressed in **nkats/µg**):

CS-11(22)	CS-11(32)	CS-11(65)
2.3 ± 0.4	3.7 ± 0.6	6.7 ± 1.0

The exact activity in nkats/µg with CS-11(22) is reported on the analysis certificate for each lot.

**Viral safety:**

The plasma used for factor Xa preparation and for BSA was tested with registered methods and found negative for bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of bovine origin must then be handled with all the required cautions, as being potentially infectious.

**Use:**

For in vitro use only. All research studies and protocols where a source of highly purified bovine Factor Xa is necessary. Studies on heparin and Low Molecular weight heparin.

**Stability:**

- Restored : - 7 days at room temperature (18-25 °C)
- >3 months at 2-8 °C
- 6 months at -20 °C or below.

Lyophilized : Until the expiration date printed on the vial.

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