Manufactured By: HYPHEN BioMed

Bovine Anti-Thrombin (AT)
Anti-Thrombine bovine (AT)
Vial of 10 mg/flacon de 10 mg

Ref. ABP104C

STORE AT
2-8 °C

For research use only

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CONSERVER A
2-8 °C

Origin:

Purified from bovine plasma using a combination of affinity chromatographies.

Presentation:

Vial containing approximately 10mg of AT corresponding to about 60 Plasma Equivalent Units or PEU (1 unit corresponds to the amount of AT in 1 ml plasma, or about 150µg), which is about 60 International Units (IU) (when determined and standardized against the WHO/NIBSC International standard for Antithrombin, plasma, human).

Reconstitution:

Each vial must be restored with 2 ml distilled water, in order to get a concentration of about 5 mg/ml, in a physiological phosphate saline buffer (pH 7.5). Dilute in the adequate buffer, if necessary.

Excipients:

Content per vial: Phosphate (16mg), Sodium Chloride (18mg).

Purity:

One major band at about 60,000 daltons on SDS-PAGE

Activity:

Tested for anti-Xa activity in presence of heparin, with a chromogenic assay (> 6IU/mg). Factor Xa neutralisation capacity is > 500 µg of Xa/mg AT.

The absence of heparin is verified after purification, using a toluidine blue assay (refer to: F.C MacIntosh, "A colorimetric method for the standardization of heparin preparations", Biochem, 1941, 35:776-782), and an anti-Xa specific chromogenic assay.

The exact activity is reported on the analysis certificate for each lot.

Viral safety:

The plasma used for AT purification was tested with registered methods and found negative for bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of bovine origin must then be handled with all the required cautions, as being potentially infectious.

Use:

For *in vitro* use only. All research studies and protocols where a source of highly purified bovine AT is necessary.

Stability:

- Restored :
- 72 hours at room temperature (18-25 °C)
 - 7 days at 2-8 °C
 - 2 months frozen at -30 °C or below

Lyophilized : Until the expiration date printed on the vial.

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