Manufactured By: HYPHEN BioMed

# Human Activated Protein C

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Vial of 100 µg/flacon de 100 µg

Ref. AEZ004C

**STORE AT**  
**2-8°C**

For research use only

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**CONSERVER A**  
**2-8°C**

**Origin:**

Highly purified activated Protein C from citrated human plasma, activated with agarose bound human thrombin.

**Presentation:**

Vial containing approximately 100 µg of highly purified human activated Protein C in presence of additives and preservatives.

**Reconstitution:**

Each vial must be restored with 1 ml distilled water. It can be diluted to the desired concentration when required.

**Excipients:**

- Glycine, hepes, Ciprofloxacin.
- 9 mg sodium chloride/vial.

**Purity:**

The Protein C used for preparing activated Protein C has one major band of 62,000 daltons on SDS-PAGE.

**Activity:**

Tested for its anticoagulant activity in an Activated Partial Thromboplastin Time assay (APTT), and for its amidolytic activity in presence or absence of PROTAC.

**Viral safety:**

The material used for Activated Protein C preparation was tested with registered methods and found negative for HIV antibodies, HBs Ag and HVC antibodies. However, no assay may warrant the total absence of infectious agents. Any product of human origin must then be handled with all the required cautions, as being potentially infectious.

**Use:**

For in vitro use only.

Activated Protein C Resistance assays.

All research studies and protocols where a source of highly purified activated Protein C is necessary.

**Stability:**

- Restored :**
- 24 hours at room temperature
  - 72 hours at 2-8 °C
  - 2 months at -20 °C or below

**Lyophilized :** - Until the expiration date printed on the vial.

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