Manufactured By: HYPHEN BioMed

Anti-FPA

Affinity purified rabbit antibodies to FPA

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CONSERVER A
2-8 °C

Origin:

Rabbit polyclonal antibodies obtained by immunisation with synthetic FPA, reproducing the sequence of human FPA (16 Amino Acids).

Antibodies are purified by immuno-affinity onto FPA coupled to agarose beads.

Presentation:

Vial containing 100 µg of rabbit antibodies specific for human FPA, affinity purified.

Reconstitution:

The vial must be reconstituted with 0.5 ml of distilled water in order to obtain an anti-FPA solution with a concentration of 200 µg/ml.

Excipients:

Glycine, Hepes, Sodium Chloride (4.5 mg per vial).

Immunogen:

Synthetic human FPA (16 Amino Acids) coupled to Bovine Serum Albumin (BSA) as carrier.

Reactivity:

Is highly reactive with human FPA, and with the FPA moieties of human fibrinogen. Does not react with BSA.

Use:

Any method where a source of highly purified anti-FPA antibodies is required.

When anti-FPA is used diluted, it is recommended to use a diluent containing 1% of BSA or of casein in order to avoid non-specific adsorption of antibodies on the recipient walls.

Stability:

Once restored, anti-FPA is stable, in its original vial, 1 month at 2-8 °C. It is recommended to add sodium azide* at 0.9 g/l, as preservative

***CAUTION:** Sodium azide (NaN₃), may react with lead and copper plumbing to form highly explosive metal azides. Flush with large volumes of water when discarding into a sink.

For prolonged conservation, it is recommended to split anti-FPA in small aliquots and to store them frozen at -20 °C, or below.

Lyophilized : Until the expiration date printed on the vial.

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