Manufactured By: HYPHEN BioMed

Purified B2-Glyco-Protein I (B2GPI)
B2-Glyco-Protéine I purifiée (B2GPI)
Vial of 100 µg/flacon de 100 µg

Ref. APP010A

STORE AT
2-8 °C

For research use only

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CONSERVER A
2-8 °C

Origin:

Highly purified human B2 Glyco Protein I (B2GPI) extracted from non activated and lipid poor human plasma, using soft and non-denaturing techniques (combination of ion-exchange chromatographies, affinity chromatographies and gel filtration).

Presentation:

Vial containing approximately 100 µg of highly purified human B2GPI, in presence of additives and preservatives.

Reconstitution:

Each vial must be restored with 1 ml distilled water. It can be diluted at the desired concentration in the appropriate buffer when required.

Nota: For using B21GPI, at low concentrations, in research studies, it is recommended to use a buffer containing bovine serum albumine or casein. This prevents protein binding to recipient walls and therefore prevents losses.

When B2GPI must be coated onto an ELISA plate, it must be diluted in a freshly prepared 0.05M Na₂HPO₄ buffer at pH 7.5. Recommended coating concentration is 10 µg/ml.

Excipients:

– Glycine, sodium chloride and hepes

Purity:

B2GPI shows a major band at 50,000 daltons when tested on SDS-PAGE.

Specific activity:

B2GPI activity is of about 4 units (one unit is the amount in 1 ml normal human plasma) per mg.

Viral safety:

The material used for B2GPI preparation was tested with registered methods and found negative for HIV antibodies, HBs Ag and HVC antibodies. However, no assay may warrant the total absence of infectious agents. Any product of human origin must then be handled with all the required cautions, as being potentially infectious.

Use:

For in vitro use only. All research studies and protocols where a source of highly purified, non degraded, human B2GPI is required. Protein used for coating microELISA plates for the assay of autoantibodies to B2GPI, present in some plasmas with anti-phospholipid antibodies.

Stability:

Restored :

- 24 hours at room temperature
- 72 hours at 2-8 °C
- 2 months frozen at -20°C or below

Lyophilized : Until the expiration date printed on the vial.

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