# BIOPHEN™ CS-41(03)

## Plasmin and Plasminogen – SK

### Chromogenic substrate

Vial of 25 mg/flacon de 25 mg

Ref. 229041

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Synthetic generic Chromogenic substrate, highly purified and stabilized.

**Specificity:**

Recommended substrate for Plasmin and Plasminogen - SK.

**Presentation:**

Vial containing 25 mg of BIOPHEN™ CS-41(03), lyophilised in presence of Mannitol as a bulking agent.

**Reconstitution:**

According to the research protocol used, the BIOPHEN™ CS-41(03) chromogenic substrate can be restored with variable volumes of distilled water; for example 5 mL can be used for a substrate concentration of 5 mg/mL, or 20 mL for a substrate concentration of 1.25 mg/mL..

**Peptide sequence:**

PyroGlu–Phe-Lys-pNa-HCl

**Purity grade:**

&gt; 95%.

**Molecular weight:**

524.6.

**pNA content (measured on RP-HPLC)**

RP-HPLC &lt; 0.1%.

**Use:**All research studies and protocols where a source of chromogenic substrate for Plasmin and Plasminogen-SK is required. **This kit is for research use only and should not be used for patient diagnosis or treatment.****Respective reactivities:**

Plasmin	Thrombin	FXa	Kallicrein	aPC
100	0	1	/	0

Assay conditions must be established for making the substrate totally specific for Plasmin and Plasminogen-SK.

**Stability:**

- Restored :**
- 7 days at room temperature
  - 3 months at 2-8 °C
  - **Do not freeze.**

**Lyophilized :** - Until the expiration date printed on the packaging.

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# BIOPHEN™ CS-41(03)

## Plasmin and Plasminogen – SK Chromogenic substrate

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DATE :

CONCLUSIONS :  Passed/Conforme  Refused/Non-conforme

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