BIOPHEN CS-01(38)

Thrombin Chromogenic substrate

Vial of 100 mg/flacon de 100 mg

Ref. 229101

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English, last revision 03-2017

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ORIGIN:

Synthetic generic Chromogenic substrate, highly purified and stabilized.

SPECIFICITY:

Recommended substrate for Thrombin.

PRESENTATION:

Vial containing about 100 mg of BIOPHEN CS-01(38), lyophilised in presence of Mannitol as a bulking agent.

This kit is for research use only and should not be used for patient diagnosis or treatment.**RECONSTITUTION:**

According to the research protocol used, the BIOPHEN CS-01(38) chromogenic substrate can be restored with variable volumes of distilled water; for example 5 mL can be used for a substrate concentration of 20 mg/mL.

PEPTIDE SEQUENCE:

D-Phe-Pip-Arg-pNa, 2HCl.

PURITY GRADE:

≥ 95%.

MOLECULAR WEIGHT:

553.

FREE PNA CONTENT (MEASURED ON SUBSTRATE AT 2.5G/L):OD_{405nm} ≤ 0.30.**USE:**

All research studies and protocols where a source of chromogenic substrate for Thrombin is required.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.**RESPECTIVE REACTIVITIES:**

Thrombin	FXa	Plasmin	Kallicrein	aPC
100	5	5	60	40

Assay conditions must be established for making the substrate totally specific for Thrombin.

STABILITY:**Restored :**

- 7 days at room temperature (18-25°C).
- 3 months at 2-8°C.
- **Do not freeze.**

Lyophilized :

- Until the expiration date printed on the vial.

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