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Français, dernière revision 06-2017

COMPOSITION:

Vial of 10 mL of Tris buffer (0.5 M) containing sodium chloride (1.75 M), Na₂ EDTA (0.075 M), PEG6000 (1%) and Sodium azide (0.9 g/L), as preservative, 10 fold concentrated (x10).

*CAUTION: Sodium azide (NaN₃), may react with lead and copper plumbing to form highly explosive metal azides. Flush with large volumes of water when discarding into a sink.

This kit is for research use only and should not be used for patient diagnosis or treatment.

PH:

8.40 (at the 1:10 dilution).

USE:

10x concentrated buffer for chromogenic assays.

Dilute at 1:10 with distilled water before use.

Once diluted, can be used as the buffer recommended by the US Pharmacopoeia for Heparin anti-Xa (eg see #221010) two-stages chromogenic assays.

Note 1: Take care of stabilizing the solution at room temperature (18-25°C) for 30 minutes before use.

Note 2: Refer to the specific associated kit inserts for instructions concerning the tests to be performed.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

STABILITY:

In its original vial, unopened and kept at 2-8°C, until the expiration date printed on the label.

When open and protected from any contamination, this solution is stable for 4 weeks at 2-8°C.

Changes compared to the previous version.

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