Recombinant Hirudin

Hirudine recombinante

Vial of $\approx 16,000$ ATU /flacon de $\approx 16\ 000$ ATU (RE120A)

RE120A

STORE AT
2-8 °C

For research use only.
Not for use in diagnostic procedures

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CONSERVER A
2-8 °C

Origin:
Recombinant protein

Presentation:
Vial containing about 16,000 ATU of rec-Hirudin.

This specification is based on the hirudin anti-thrombin activity expressed in ATU*.

Reconstitution:
Restore the vial with 1 ml of distilled water. When restored, the solution contains a rec-Hirudin solution with an activity of approximately 16,000 ATU*/ml.
The measured recombinant hirudin activity is given for each lot on the Certificated of Analysis.

Excipients:
– Glycine, Hepes , Sodium Chloride, **pH 7.50**

Activity:
Hirudin has an activity of about 16,000 ATU*/vial

Use:
For in vitro use only.
For research use only (RUO)
Thrombin inhibitor, which can be used in any in vitro laboratory research application where an efficient inhibition of Thrombin is required.

Stability:
Restored:
– 2 days at room temperature
– 7 days at 2-8 °C
– 6 months frozen at -20 °C or below

Lyophilized :
– Until the expiration date printed on the vial.

*ATU : Anti Thrombin Unit

This kit is for research use only and should not be used for patient diagnosis or treatment. The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

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