Recombinant human tPA

tPA humain recombinant

Vial of 10 µg/flacon de 10 µg

Ref. RP007A

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Recombinant human tPA.

PRESENTATION:

Vial containing 10 µg of recombinant human tPA, lyophilized.

This kit is for research use only and should not be used for patient diagnosis or treatment.**RECONSTITUTION:**

Restore the vial with 1 mL of distilled water.

Nota : If tPA must be used at a concentration below 0.01 mg/mL, the dilution buffer must contain Bovine Serum Albumin (i.e. 1%) or any other carrier protein or polymer in order to avoid loss of the protein by adsorption on the container walls.

EXCIPIENTS:Phosphate, Arginine, Sodium Chloride, pH 7.50.
Bovine Serum Albumin 1%.**VIRAL SAFETY:**

The bovine plasma used for BSA preparation was tested with registered methods and found negative for bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of biological origin must then be handled with all the required cautions, as being potentially infectious.

ACTIVITY:

Recombinant human tPA has an activity of about 500,000 IU/mg (tested for its capacity to induce a fibrin clot lysis). tPA activity is labile. Recombinant human tPA must be used rapidly or stored frozen in small aliquots and thawed just before use.

CHARACTERISTICS:Recombinant human tPA has a molecular weight of about 59,000 daltons.
Recombinant human tPA preparation has a purity grade > 95%.**USE:**For *in vitro* use only.

All research studies and protocols where a source of recombinant human tPA is necessary.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.**STABILITY:**

Restored :

- 12 hours at room temperature (18-25°C).
- 24 hours at 2-8°C.
- 6 months at -20°C or below.

Lyophilized :

- Until the expiration date printed on the vial.

Freezing conditions and stability of the thawed product should be checked in the working conditions of the laboratory user.

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