

ANIARA

Manufactured By: Medicago AB

Antiserum against *Encephalitozoon cuniculi*



Features

- Antiserum against *E. cuniculi*
- Different animal sources available (guinea pig, rabbit, rat)

Product description

The serum is produced from naturally-infected animals. 0.05% azide is added. Before use in a test, dilute the serum 1:20 or more in PBS.

For laboratory use only.

Applications

- Veterinary procedures
- A parasite reagent that can be used as a positive control in immunoassays

Directions for use

Dilute serum 1:20 or more in PBS buffer. Handle as non-infectious. Use lab coat, gloves and goggles.

Shipping and storage

The product is shipped at room temperature, store diluted serum at 2°C to 8°C.

Certifications

Medicago's laboratories and manufacturing site in Uppsala are ISO 9001:2008 and ISO 13485:2003 certified. Each stage of the manufacturing process is controlled and monitored by stringent quality control procedures to guarantee the highest possible quality and lot-to-lot reproducibility.



Specifications

Content	Serum from Rabbit/Rat/Guinea pig immunized with <i>E.cuniculi</i>
Formulation	0.05% azide (NaN ₃)
Storage	2°C to 8°C
Shelf life	12 months after dilution

Ordering information

Article no.	Product name	Pack size
A18-4004-1	Guinea pig antiserum against <i>E. cuniculi</i>	1 ml
A18-4001-1	Rabbit antiserum against <i>E.cuniculi</i>	1 ml
A18-4002-1	Rat antiserum against <i>E.cuniculi</i>	1 ml

v. 01 2010-08