# BIOPHEN CS-11(22)

## Factor Xa Chromogenic Substrate

Vial of 25 mg/flacon de 25 mg

Ref. 229015

**STORE AT**  
**2-8 °C**

**For research use only**  
**Not for use in diagnosis**  
**procedures.**

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**CONSERVER A**  
**2-8 °C**

**Origin:**

Synthetic generic Chromogenic substrate, highly purified and stabilized.

**Specificity:**

Recommended substrate for Factor Xa.

**Presentation:**

Vial containing 25 mg of BIOPHEN CS-11(22), lyophilised in presence of Mannitol as a bulking agent.

**Reconstitution:**

According to the research protocol used, the BIOPHEN CS-11(22) chromogenic substrate can be restored with variable volumes of distilled water ; for example 10 mL can be used for a substrate concentration of 2.5 mg/mL, or 20 mL for a substrate concentration of 1.25 mg/mL. Shake thoroughly until complete dissolution (vortex). Let to stabilize for 30 min. at room temperature.

**Peptide sequence:**

Mixture (50%-50%) of Bz-Ile-Glu ( $\gamma$ OCH<sub>3</sub>)-Gly-Arg-pNa (form 1) and Bz-Ile-Glu ( $\gamma$ OH)-Gly-Arg-pNa (form 2)

**Purity grade:**

> 95%.

**Molecular weight:**

711.8 (form 1) and 697.7 (form 2).

**Free pNA content (measured on substrate at 2.5g/L):**

OD<sub>405nm</sub> ≤ 0.30

**Use:**

All research studies and protocols where a source of chromogenic substrate for Factor Xa is required. **This kit is for research use only and should not be used for patient diagnosis or treatment.**

**Respective reactivities:**

FXa	Thrombin	Plasmin	Kallicrein	aPC
100	1	2	3	0

Assay conditions must be duly established for rendering the assay conditions totally specific for Factor Xa, when this substrate is used.

**Stability:**

**Restored :**

- 7 days at room temperature
- 3 months at 2-8 °C
- **Do not freeze.**

**Lyophilized :** Until the expiration date printed on the vial.

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