Purified bovine Factor Xa

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Vial of 30 µg or set of 10x30 µg
flacon de 30 µg ou coffret de 10x30 µg

Ref. BE101D/BE101L

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English, last revision 12-2016

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ORIGIN:

Highly purified Factor X from citrated bovine plasma, activated with agarose bound RVV.

PRESENTATION:

Each Vial containing approximately 30 µg of highly purified bovine Factor Xa in presence of additives and preservatives.

(1 Unit FXa is the amount which can be generated from 1 mL citrated bovine plasma, ie about 10 µg).

This kit is for research use only and should not be used for patient diagnosis or treatment.

RECONSTITUTION:

Each vial must be restored with 10 mL distilled water, in order to get a concentration of about 3 µg/mL. It can be diluted to the desired concentration, when required, in the adequate buffer.

EXCIPIENTS:

Content per vial: Tris (60mg) , Sodium chlorure (50mg), PEG 6000 (10mg), BSA (20mg).

PURITY:

The Factor X zymogen used for preparing Factor Xa has one major band of about 55,000 daltons on SDS-PAGE.

ACTIVITY:

Tested for its amidolytic activity on a FXa specific substrate, in a purified system without or with RVV. All FX is converted to FXa.

Chromogenic activity on specific substrates (expressed in nkats): tested in the optimized conditions (0.05M Tris buffer at pH8.40, containing 0.30M NaCl), the specific activities are defined on these FXa substrates:

CS-11(22)	CS-11(32)	CS-11(65)
-----------	-----------	-----------

The exact activity in nkats/µg is reported on the analysis certificate for each substrate.

VIRAL SAFETY:

The plasma used for Factor Xa preparation and for BSA was tested with registered methods and found negative for bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of bovine origin must then be handled with all the required cautions, as being potentially infectious.

USE:

For *in vitro* use only. All research studies and protocols where a source of highly purified bovine Factor Xa is necessary. Studies on heparin and Low Molecular weight heparin.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

STABILITY:

Restored : - 7 days at room temperature (18-25°C).
- 3 months at 2-8 °C.
- 6 months at -20°C or below.
Lyophilized : Until the expiration date printed on the kit.

The stability of the reconstituted reagent should be checked under laboratory work conditions.

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