Purified bovine Factor Xa

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Vial of 50 µg or set of 6x50 µg
flacon de 50 µg ou coffret de 6x50 µg

Ref.BE1010/BE101K

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English, last revision 12-2016

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ORIGIN:

Highly purified Factor X from citrated bovine plasma, activated with agarose bound RVV.

PRESENTATION:

Each vial containing approximately 50 µg of highly purified bovine Factor Xa in presence of additives and preservatives.
(1 Unit FXa is the amount which can be generated from 1 mL citrated bovine plasma, ie about 10 µg).
This kit is for research use only and should not be used for patient diagnosis or treatment.

RECONSTITUTION:

Each vial must be restored with 1 mL distilled water, in order to get a concentration of about 50µg/mL. It can be diluted to the desired concentration, when required, in the adequate buffer.

EXCIPIENTS:

Content per vial: Glycine (20mg), Hepes (6mg), Prionex (0.25% after reconstitution), BSA (10mg), Sodium chloride (9mg).

PURITY:

The Factor X zymogen used for preparing Factor Xa has one major band of about 55,000 daltons on SDS-PAGE.

ACTIVITY:

- a) Tested for its amidolytic activity on a FXa specific substrate, in a purified system without or with RVV. All FX is converted to FXa.
b) Chromogenic activity on specific substrates (expressed in nkats): tested in the optimized conditions (0.05M Tris buffer at pH8.40, containing 0.30M NaCl), the specific activities are defined on these FXa substrates:

CS-11(22)	CS-11(32)	CS-11(65)
-----------	-----------	-----------

The exact activity in nkats/µg is reported on the analysis certificate for each substrate.

VIRAL SAFETY:

The plasma used for Factor Xa preparation and for BSA was tested with registered methods and found negative for bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of bovine origin must then be handled with all the required cautions, as being potentially infectious.

USE:

For *in vitro* use only. All research studies and protocols where a source of highly purified bovine Factor Xa is necessary. Studies on heparin and Low Molecular weight heparin.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

STABILITY:

Restored : - 30 days at room temperature (18-25°C).
- 3 months at 2-8 °C.
- 6 months at -20°C or below.
Lyophilized : Until the expiration date printed on the vial.

The stability of the reconstituted reagent should be checked under laboratory work conditions.

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