Human Activated Protein C

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Vial of 10 µg/flacon de 10 µg

Ref. EZ004B

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English, last revision 03-2017

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ORIGIN:

Highly purified activated Protein C from citrated human plasma, activated with agarose bound human thrombin.

PRESENTATION:

Vial containing approximately 10 µg of highly purified human activated Protein C in presence of additives and preservatives.

This kit is for research use only and should not be used for patient diagnosis or treatment.

RECONSTITUTION:

Each vial must be restored with 1 mL distilled water. It can be diluted to the desired concentration when required.

EXCIPIENTS:

Tris, PEG 6000, Bovine Albumin Serum, Ciprofloxacin.
9 mg sodium chloride/vial.

PURITY:

The Protein C used for preparing activated Protein C has one major band of 61,000 daltons on SDS-PAGE.

ACTIVITY:

Tested for its anticoagulant activity in an Activated Partial Thromboplastin Time assay (aPTT), and for its amidolytic activity in presence or absence of PROTAC.

VIRAL SAFETY:

The material used for Activated Protein C preparation was tested with registered methods and found negative for HIV antibodies, HBs Ag, HCV antibodies and bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of human or bovine origin must then be handled with all the required cautions, as being potentially infectious.

USE:

For *in vitro* use only.

Activated Protein C Resistance assays.

All research studies and protocols where a source of highly purified activated Protein C is necessary.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

STABILITY:

Restored :

- 24 hours at room temperature (18-25°C).
- 72 hours at 2-8°C.
- 2 months at -20°C or below.

Lyophilized :

- Until the expiration date printed on the vial.

Freezing conditions and stability of the thawed product should be checked in the working conditions of the laboratory user.

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