Purified human Factor Xa

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Vial of 10 µg / Flacon de 10 µg
Set of 6x10 µg / Coffret de 6x10 µg

Ref. EZ007A / EZ007K

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English, last revision: 12-2016

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ORIGIN:

Highly purified Factor X from citrated human plasma, activated with agarose bound RVV.
MW of about 44 000.

PRESENTATION:

Vial containing approximately 10 µg of highly purified human Factor Xa in presence of additives and preservatives.
(1 Unit FXa is the amount which can be generated from 1 mL citrated human plasma, ie about 10 µg).

This kit is for research use only and should not be used for patient diagnosis or treatment.

RECONSTITUTION:

Each vial must be restored with 1 mL distilled water, in order to get a concentration of about 10µg/mL. It can be diluted to the desired concentration, when required, in the adequate buffer.

EXCIPIENTS:

Content per vial: Glycine (20mg), Hepes (6mg), Prionex (0.5% after reconstitution), PEG6000 (5 mg), Sodium chloride (9mg).

PURITY:

The Factor X used for preparing Factor Xa has one major band of 59,000 daltons on SDS-PAGE.

ACTIVITY:

a) Tested for its amidolytic activity on a FXa specific substrate, in a purified system without or with RVV. All FX is converted to FXa.
b) Chromogenic activity on specific substrates (expressed in nkats) tested in the optimized conditions (0.05M Tris buffer at pH8.40, containing 0.30M NaCl), the specific activities are defined on various FXa substrates :

CS-11(22)	CS-11(32)	CS-11(65)
-----------	-----------	-----------

The exact activity in nkats/µg with CS-11(32) is reported on the analysis certificate for each lot.

VIRAL SAFETY:

The human plasma used for Factor Xa preparation was tested with registered methods and found negative for HIV (1 and 2) antibodies, HBs Ag and HVC antibodies. However, no assay may warrant the total absence of infectious agents. Any product of human origin must then be handled with all the required cautions, as being potentially infectious.

USE:

For in vitro use only. All research studies and protocols where a source of highly purified human Factor Xa is necessary. Studies on heparin and Low Molecular weight heparin.
The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

STABILITY:

Restored:

- 7 days at room temperature (18-25°C).
- 1 month at 2-8 °C.
- 6 months at -20°C or below.

Lyophilized: Until the expiration date printed on the vial.

The stability of the reconstituted reagent should be checked under laboratory work conditions.

Changes compared to the previous version.

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