**Protac<sup>®</sup>**

Vial of 5 units/flacon de 5 unités

Ref. EZ011B

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English, last revision 03-2017

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**ORIGIN:**Protac<sup>®</sup>, Protein C Activator extracted from the snake venom of Agkistrodon contortrix, purified and stabilised.**PRESENTATION:**Vial containing approximately about 5 Units of Protac<sup>®</sup> lyophilised in presence of additives and preservatives.**This kit is for research use only and should not be used for patient diagnosis or treatment.****RECONSTITUTION:**

Each vial must be restored with 1 mL distilled water in order to obtain a preparation at 5 U/mL.

**EXCIPIENTS:**

Glycine, Hepes, Ciprofloxacin and Bovine Serum Albumin (BSA).

**PURITY:**Protac<sup>®</sup> is a highly purified enzyme from Agkistrodon contortrix, with a major band of 36 to 42 Kdaltons on SDS-PAGE.**ACTIVITY:**Protac<sup>®</sup> is tested for its capacity to activate the Protein C into Activated Protein C (APC). Tested with chrometric and chromogenic methods.**VIRAL SAFETY:**Protac<sup>®</sup> preparation contains BSA, which was tested with registered methods and found negative for bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of bovine origin must then be handled with all the required cautions, as being potentially infectious.**Cautions: Protac<sup>®</sup> is extracted from snake venom. Avoid contact with skin and eyes. In case of contact flush with copious amount of water. Handle this product with all the required cautions for this kind of reagents.****USE:**For *in vitro* research use only.**The results obtained should be for research purposes only and not used for patient diagnosis or treatment.****STABILITY:**

Restored :

- 24 hours at room temperature (18-25°C).
- 72 hours at 2-8°C.
- 2 months at -20°C or below.

Lyophilized :

- Until the expiration date printed on the vial.

Freezing conditions and stability of the thawed product should be checked in the working conditions of the laboratory user.

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DATE :

CONCLUSIONS :  Passed/Conforme  Refused/Non-conforme

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