# Purified human plasmin

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Vial of 2.5 IU / Flacon de 2,5 UI

Ref. EZ014A

**STORE AT**  
**2-8 °C**

**FOR RESEARCH USE ONLY.**  
**NOT FOR USE IN DIAGNOSTIC**  
**PROCEDURES.**

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English, last revision 03-2017

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**ORIGIN:**

Purified plasmin from citrated human plasma.

**PRESENTATION:**

Vial containing approximately 2.5 IU of purified human plasmin in presence of additives and preservatives.

**This kit is for research use only and should not be used for patient diagnosis or treatment.**

**RECONSTITUTION:**

Vial must be restored with 1 mL distilled water. It can be diluted to the desired concentration, when required, in the adequate buffer. For a better stability, vials can be restored with 1 mL of distilled water containing 25% glycerol.

**EXCIPIENTS:**

Bovine Albumin Serum (1%), Mannitol, preservatives.

**PURITY:**

The plasminogen used for preparing plasmin has one major band of 88,000 daltons on SDS-PAGE.

**ACTIVITY:**

2.5 IU.

**VIRAL SAFETY:**

The human plasma used for human plasmin preparation was tested with registered methods and found negative for HIV (1 and 2) antibodies, HBs Ag, HCV antibodies and bovine infectious diseases, notably for the bovine spongiform encephalopathy. However, no assay may warrant the total absence of infectious agents. Any product of human or bovine origin must then be handled with all the required cautions, as being potentially infectious.

**USE:**

This product is for research use only and should not be used for patient diagnosis or treatment. All research studies and protocols where a source of highly purified human plasmin is necessary.

**The results obtained should be for research purposes only and not used for patient diagnosis or treatment.**

**STABILITY:**

Lyophilized:

- Until the expiration date printed on the vial.

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