# Human Protein S

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Vial of 100 µg/flacon de 100 µg

Ref. PP012A

**STORE AT**  
**2-8 °C**

For research use only

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**CONSERVER A**  
**2-8 °C****Origin:**

Highly purified Protein S extracted from citrated human plasma.

**Presentation:**

Vial containing approximately 100 µg of highly purified human Protein S in presence of additives and preservatives.

**Reconstitution:**

Each vial must be restored with 1 mL distilled water.

**Excipients:**

– Glycine, sodium chloride, hepes.

**Purity:**

One major band of 69,000 daltons on SDS-PAGE

**Specific activity:**

Tested on SDS-PAGE, for its anticoagulant activity as activated Protein C cofactor, for the antigen concentration and for the absence of factor II in a factor II deficient plasma system.

≥ 8 PEU/vial of Protein S activity (1 Plasma Equivalent Unit or 1 PEU is the activity in 1 mL of normal human plasma)

**Viral safety:**

The plasma used for Protein S purification was tested with registered methods and found negative for HIV antibodies, HBs Ag and HVC antibodies. However, no assay may warrant the total absence of infectious agents. Any product of human origin must then be handled with all the required cautions, as being potentially infectious.

**Use:**For *in vitro* use only.

All research studies and protocols where a source of highly purified Protein S is necessary.

**Stability:****Restored :**  
- 72 hours at room temperature (18-25 °C)  
- 3 days at 2-8 °C  
- 6 months frozen at -20 °C or below**Lyophilized :** Until the expiration date printed on the vial.**Origine :**

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CONSERVER A  
2-8 °C

## ANALYSIS CERTIFICATE

ANALYTICAL DATA	SPECIFICATIONS
<b>1. Protein Concentration (Lowry method)</b>  <div style="text-align: right;">..... µg/vial</div>	>90 µg/vial
<b>2. Protein S antigenic concentration</b>  <b>Free</b> ..... U/vial                      ..... U/mg <b>Total</b> ..... U/vial                      ..... U/mg	≥ 8 U/vial ( <b>Free</b> ) ≥ 3 U/vial ( <b>Total</b> )
<b>3. SDS-PAGE (5 % acrylamide)</b>  <div style="text-align: center;">Protein S has a major band of about ..... daltons</div>	1 major band of about 69,000 daltons
<b>4. Protein S activity</b>  <div style="text-align: right;">..... U/vial</div>	≥6 U/vial
<b>5. Prothrombin concentration</b>  <div style="text-align: right;">&lt; ...../vial</div>	≤ 0.01 U/vial

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DATE :

CONCLUSIONS :  Passed/Conforme  Refused/Non-conforme

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