# Heparin cofactor II

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Vial of 1 mg/flacon de 1 mg

Ref. PP014C

**STORE AT**  
**2-8 °C**

For research use only

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**CONSERVER A**  
**2-8 °C**

**Origin:**

Highly purified Heparin Cofactor II extracted from human plasma using a combination of salt and PEG precipitation and affinity and ion-exchange chromatographies.

**Presentation:**

Vial containing approximately 1 mg of heparin Cofactor II corresponding to about 10 Plasma Equivalent Units or PEU (amount in 1 mL human plasma).

**Reconstitution:**

Each vial must be restored with 1 mL distilled water.

**Excipients:**

- Phosphate buffer
- 9 mg sodium chloride/vial

**Purity:**

One major band at 66,000 daltons on SDS-PAGE

**Activity:**

Tested in a chromogenic assay for the thrombin inhibition capacity in presence of an excess of dermatan sulphate.

**Viral safety:**

The plasma used for Heparin Cofactor II purification was tested with registered methods and found negative for HIV antibodies, HBs Ag and HVC antibodies. However, no assay may warrant the total absence of infectious agents. Any product of human origin must then be handled with all the required cautions, as being potentially infectious.

**Use:**

For *in vitro* use only.  
All research studies and protocols where a source of highly purified Heparin Cofactor III is necessary.

**Stability:**

- Restored :**
- 24 hours at room temperature (18-25 °C)
  - 72 hours at 2-8 °C
  - 2 months frozen at -20 °C or below

**Lyophilized :** Until the expiration date printed on the vial.

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