Rabbit Recombinant Tissue Factor

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English, last revision 03-2017

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PRESENTATION:

Vial containing approximately 50 µg of Rabbit recombinant Tissue Factor lyophilised in presence of additives and preservatives.

This kit is for research use only and should not be used for patient diagnosis or treatment.

RECONSTITUTION:

Each vial must be restored with 1 mL distilled water.

EXCIPIENTS:

Tris 0.05M, NaCl 0.15M, Mannitol 2%, CHAPS 0.01%, pH 7.50.

PURITY:

One major band of about 46,000 daltons on SDS-PAGE.

SPECIFIC ACTIVITY:

Tested for its clotting activity, in activating FVII in presence of calcium, when relipidated.

USE:

For *in vitro* use only.

All research studies and protocols where a source of Rabbit recombinant Tissue Factor is necessary.

Tissue Factor clotting activity is fully expressed when it is appropriately relipidated, and tested in presence of calcium.

The results obtained should be for research purposes only and not used for patient diagnosis or treatment.

STABILITY:

Restored :

- 24 hours at room temperature (18-25°C).
- 3 days at 2-8°C.
- 6 months frozen at -20°C or below.

Lyophilized :

- Until the expiration date printed on the vial.

Freezing conditions and stability of the thawed product should be checked in the working conditions of the laboratory user.

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