



Manufactured By: Medicago AB

Immunosorb A



Features

- Pure recombinant Protein A
- Immobilized on a cross-linked agarose matrix
- Strong binding to Fc region of IgG
- High capacity: > 20 mg IgG/ml

Product description

Immunosorb A consists of pure recombinant protein A which has been derived from the bacterium *Staphylococcus aureus*. The protein is cloned and expressed in *E.coli* and immobilized on a cross-linked agarose matrix.

Protein A binds to specific classes of immunoglobulins and can preferably be used for purification of mouse and rabbit immunoglobulins. The protein binds to the Fc region of immunoglobulins by interacting with the heavy chain. It binds with high affinity to human IgG₁ and IgG₂ as well as mouse IgG_{2a} and IgG_{2b}. Protein A binds to human IgM, IgA and IgE as well as to mouse IgG₃ and IgG₁ but with moderate affinity. It does not react with human IgG₃ or IgD, nor with mouse IgM, IgA or IgE.

The material is non-cultural, non-infectious and is not pathogenic for animals or plants. It does not contain any animal or cell culture derived products or additions such as albumin or serum nor any toxic or harmful substances.

Immunosorb A is provided as a liquid solution with 20 % ethanol as preservative.

For research use only.

Applications

- Purification of mouse and rabbit immunoglobulins
- Immobilization and detection of immunoglobulins
- Affinity chromatography

Specifications

Content	Recombinant Protein A
Format	White liquid solution, 20 % EtOH
Source	<i>Staphylococcus aureus</i>
Purity	> 95 %
Package	flasks
Storage	4°C

Directions for use

No additional mixing reagents are required. The material is not dependent on room temperature.

Shipping and storage

The product is shipped at room temperature and is stable for 12 months if stored at 2°C to 8°C.

Ordering information

Article no.	Product name	Pack size
A10-1257-200	Immunosorb A	200 ml
A10-1257-100	Immunosorb A	100 ml
A10-1257-5	Immunosorb A	5 ml



7768 Service Center Drive • West Chester OH 45069

Phone: 513.770.1991

Toll Free: 866.783.3797

Fax: 513.573.9241

Email: info@aniara.com

www.aniara.com