

**U-CyTech BV**

Yalelaan 48  
3584 CM Utrecht  
The Netherlands  
P +31.30.253 5960  
F +31.30.253 9344  
INFO@ucytech.com  
www.ucytech.com

## Data sheet Monkey Perforin ELISPOT antibody pair; 20-plate format

Cat. No.: CT620-20

### Coating antibodies (1 vial)

Product: Monoclonal antibody to monkey perforin  
Isotype: Mouse IgG<sub>1</sub>  
Production: *In vitro* using serum free medium  
Purification: DEAE ion exchange chromatography  
Contents: Each vial contains sufficient material for coating of twenty 96-well ELISPOT plates  
Buffer: Prior to lyophilization: 1.0 ml PBS + 125 mM trehalose  
Application: Coating antibody in an ELISPOT system  
Reconstitution: Dissolve the contents of the vial by injection of 1.0 ml distilled water into the vial and dilute 100 times in PBS. The total amount of one vial is sufficient for twenty 96-well ELISPOT plates (1920 determinations; 50 µl/well).

### Detection antibodies (1 vial)

Product: Biotinylated monoclonal antibody to monkey perforin  
Isotype: Mouse IgG<sub>1</sub>  
Production: *In vitro* using serum free medium  
Purification: DEAE ion exchange chromatography  
Labeling: With Biotin-7-NHS (N-hydroxysuccinimide)  
Contents: Each vial contains sufficient material for twenty 96-well ELISPOT plates  
Buffer: Prior to lyophilization: 2.0 ml PBS + 1% BSA + 125 mM trehalose  
Application: Detection antibody in an ELISPOT system  
Reconstitution: Dissolve the contents of the vial by injection of 2.0 ml distilled water into the vial and dilute 100 times in Dilution buffer (see Technical Data Sheet). The total amount of one vial is sufficient for twenty 96-well ELISPOT plates (1920 determinations; 100 µl/well).

### General

Specificity: Validated for detecting rhesus macaque and barbary macaque perforin  
Sterility: Membrane filtered (0.2 µm)  
Stability: The lyophilized products are stable for at least one year at 4°C (expiry date is indicated on the vials).  
After reconstitution, the antibodies are stable for several months at 4°C (if kept sterile) or for minimal one year at -20°C.

PROUDLY DISTRIBUTED BY

**ANIARA**  
ANIARA DIAGNOSTICA LLC

+1 (513) 770-1991 | +1 (866) 783-3797  
7768 Service Center Drive  
West Chester, OH 45069  
info@aniara.com