

## Data sheet Human Perforin ELISPOT antibody pair; 20-plate format

Cat. No.: CT653-20

### Coating antibodies (1 vial)

Product: Monoclonal antibody to human perforin  
Isotype: Mouse IgG<sub>1</sub>  
Production: *In vitro* using serum free medium  
Purification: DEAE ion exchange chromatography  
Contents: Each vial contains sufficient material for coating of twenty 96-well ELISPOT plates  
Buffer: Prior to lyophilization: 1.0 ml PBS + 125 mM trehalose  
Application: Coating antibody in an ELISPOT system  
Reconstitution: Dissolve the contents of the vial by injection of 1.0 ml distilled water into the vial and dilute 100 times in PBS. The total amount of one vial is sufficient for twenty 96-well ELISPOT plates (1920 determinations; 50 µl/well).

### Detection antibodies (1 vial)

Product: Biotinylated monoclonal antibody to human perforin  
Isotype: Mouse IgG<sub>1</sub>  
Production: *In vitro* using serum free medium  
Purification: DEAE ion exchange chromatography  
Labeling: With Biotin-7-NHS (N-hydroxysuccinimide)  
Contents: Each vial contains sufficient material for twenty 96-well ELISPOT plates  
Buffer: Prior to lyophilization: 2.0 ml PBS + 1% BSA + 125 mM trehalose  
Application: Detection antibody in an ELISPOT system  
Reconstitution: Dissolve the contents of the vial by injection of 2.0 ml distilled water into the vial and dilute 100 times in Dilution buffer (see Technical Data Sheet). The total amount of one vial is sufficient for twenty 96-well ELISPOT plates (1920 determinations; 100 µl/well).

### General

Specificity: Validated for detecting human perforin  
Sterility: Membrane filtered (0.2 µm)  
Stability: The lyophilized products are stable for at least one year at 4°C (expiry date is indicated on the vials).  
After reconstitution, the antibodies are stable for several months at 4°C (if kept sterile) or for minimal one year at -20°C.

For research use only

PROUDLY DISTRIBUTED BY   
ANIARA DIAGNOSTICA LLC  
+1 (513) 770-1991 | +1 (866) 783-3797  
7768 Service Center Drive  
West Chester, OH 45069  
info@aniara.com

Version 111031