

U-CyTech BV

Yalelaan 48
3584 CM Utrecht
The Netherlands
P +31.30.253 5960
F +31.30.253 9344
INFO@ucytech.com
www.ucytech.com

Data sheet Blocking stock solution (10x)**Cat No.:** **CT363**

- Specificity:** This solution has been developed to prevent non-specific binding of detector antibodies and conjugate in the ELISPOT. The solution is suited for both the enzymatic staining procedure on PVDF membrane-bottomed plates and the silver staining procedure on polystyrene-bottomed plates.
- Application:** Prior to use, dilute the solution 10 times with pyrogen-free PBS. Apply 200 µl/well under aseptic conditions. Cover the plate and incubate for 2 hours at room temperature.
- Contents:** 10 ml.
Preservative free.
- Sterility:** Membrane filtered (0.2 µm).
- Packaging:** 10 ml clear glass bottle.
- Storage:** Refrigerated.
The solution can be safely stored at 4°C for a defined length of time (expiry date indicated on vial). After opening, the product expires after 6 months when kept sterile.
- General:** A clear light yellow solution.
Do not use if the solution is or has become turbid.
- Safety:** The product is not classified as hazardous according to Regulation (EC) no. 1272/2008 and its amendments.
Consult the material safety data sheet (MSDS) for further information regarding hazards and safe handling practices. The MSDS can be found on our website www.ucytech.com.

